

The Nelldorn Glassware Dryer



The Nelldorn Dryer provides an economical and efficient method of drying and storing most containers, glassware and plasticware. In the laboratory, volumetric flasks, measuring cylinders, test tubes and similar equipment can conveniently be dried and then left on the tubes until required. For pharmacies, the unit can be used for most types of dispensing containers including plastic.

- Cost efficient
- Capacity of 60 items
- Dries in 15 to 20 minutes
- Wall mounting or bench standing
- Thermostatic control
- Variable 120 minute timer
- Quick and Efficient

The unit will dry up to 60 items at one time. The average drying time is 15 to 20 minutes although this varies from load to load. A large volume of air is drawn over the thermostatically controlled heating elements and forced through each drying tube. This rapid flow of warm air dries glass or plastic quickly without harm. The process uses a minimum of electricity (typically less than 2 units) and is therefore considerably cheaper than oven type dryers.

Space saving design

The stepped design allows the removal of any item without disturbing the remainder of the load. This also facilitates the run-off of water to the front of the unit where a small drain type prevents it collecting the base. Air intake is recessed at the back of the unit to allow siting on any bench or even mounting on a wall.

Economic Operation

The dryer is fitted with 60 stain resistant nylon drying tubes which are rigid enough to hold position yet resilient to avoid damage to items being dried. The unit has an adjustable timer which will automatically switch off the dryer after the pre-set time from 0-120 minutes. Rocker switches control two independent 1kw cut-outs. The tangential fan blower unit is designed to give quiet, trouble free running.

DRYING TUBE CAPACITY

75mm	100mm	140mm	152mm
30	13	9	8

Total Number of Tubes	60
Voltage	220/240V, 13A
Dimensions (wxdxh)	490 X 190 430mm

For more information on The Nelldorn Dryer, contact:

Astell Scientific Ltd.

Powerscroft Road, Sidcup, Kent. DA14 5DT. UK

Tel. +44 (0) 208 309 2024/3/2 Email info@astell.com www.astell.com



Model Ref. DCU012